

THROW AWAY YOUR TIMERS!

NEVER BABYSIT YOUR SAMPLES AGAIN

THE FREEDOM ROCKER™



Model DX4 can process Westerns in up to 4 trays from blocking to final rinse.

The FREEDOM ROCKER™ automatically process blots and gels, including Westerns from blocker to final rinse. It is a rocker with a built-in pump and timer for transferring liquids to and from the trays. It frees you up from the tedious task of babysitting the blots or gels.

AUTOMATED PROCESSING ...

- > INCREASE PRODUCTIVITY
- > REDUCE DOWNTIME
- > SAVE MONEY

> INCREASE PRODUCTIVITY:

Make better use of your time. THE FREEDOM ROCKER™ automates the processing of blots and gels, freeing up time for more demanding activities. With accurate timing and pumping, results are more consistent.

> REDUCE DOWNTIME AND SHORTEN CYCLES:

THE FREEDOM ROCKER™ allows you to process blots and gels anytime ... even during meetings and over weekends.

> SAVE MONEY: Through increased productivity, THE FREEDOM ROCKER™ saves you money every time you use it. It quickly pays for itself in just 30 uses.

RISK FREE: THE FREEDOM ROCKER™ comes with a 30 day money back guarantee

"I'm one of the first users of THE FREEDOM ROCKER™ at Amgen in Cambridge, MA. The machine gets LOTS of use from all of us here." — *Dr. Barbara Grubinska, Amgen Inc.*

"It is perfect for silver staining of gels for mass spectrometry. It has made an UNBELIEVABLE impact on data generation for me in cancer protein screening." — *Dr. Steve Griffiths, Atlantic Cancer Research Institute*

"THE FREEDOM ROCKER™ is really convenient. I used it for washing gels while I worked on other things, and also as a regular rocker." — *Dr. Michele Lastro, The UAlbany Gen*NY*Sys Center for Excellence in Cancer Genomics*

NEXT>>>ADVANCE

BETTER WORKFLOW • ACCURATE RESULTS • HIGH VALUE
Next Advance, Inc. www.nextadvance.com

24 Prospect Avenue
Averill Park, NY 12018, USA
Telephone: 1.800.738.1681
Fax: 1.610.862.1718

THROW AWAY YOUR TIMERS!

NEVER BABYSIT YOUR SAMPLES AGAIN

AUTOMATICALLY PROCESS GELS/BLOTS.

The FREEDOM ROCKER™DX automates the processing of gels and blots, using your same current protocols. Programming is extremely intuitive, with each program step the same as each protocol step. Program up to 9 steps, where each step is an immersion in a reagent or a series of washes. For each step, simply set the volume and duration of the reagent or wash buffer, the number of dispenses, the rocking speed, and whether you want a clearing rinse. After the final wash, the buffer remains in the trays so that the gel or blot does not dry out.

The FREEDOM ROCKER™DX can even be programmed for non-traditional protocols requiring initial rinses. It has the capacity to fill trays from 3 sources (one at a time to all trays), and to individually recover 2 reagents as well as aspirate into a waste container or sink.

Rocking speed up to 40 cycles per minute, adjustable for each individual step, reagent and buffer immersion times from 1 minute up to 90 hours, volumes from under 10 mL to up to 600 mL.

Features:

- Unattended dispensing and aspirating of liquids.
- Can recover reagents.
- Occupies only 10 in. or 12 in. of lab bench by 14 in. deep. Platforms are 10 inch wide x 12 inch deep [25 x 30 cm, model DX2 for 1 or 2 trays] or 12 x 12 inch [30 x 30 cm, model DX4 for 1, 2, or 4 trays]. Single tray - maximum size is as large as the platform. Four trays - maximum size is large enough for minigels.
- Can use anywhere. Quiet. No vacuum or air connections or vacuum pumps required. Just plug into an electrical outlet.
- Optional clearing rinse between washes.
- Drip proof quick-release connections.
- Patent pending technology.
- CE/UL certified worldwide compatible 12V power supply.
- Customization available (e.g. alternate bottles).
- 2 year warranty.
- 30 day money back guarantee.

Model	Platform	Capacity
DX2	10" wide by 12" deep	1 or 2 Trays
DX4	12" wide by 12" deep	4 Trays; 1 or 2 Trays w/ accessory kit



BULLET BLENDER™

Lyse/Disrupt up to 24 samples at once, in just minutes. No worries about cross-contamination, sample loss, or sample heating.

OTHER INNOVATIVE PRODUCTS



SYRINGE PUMPS

Very easy to use and program. Loaded with features. Versions for up to 1, 6, or 8 syringes.



CAPILLARY LOADER

Pack HPLC capillaries. Load Mass Spec samples. Available with integrated magnetic stirplate.



NEXT ROCKERS

Very quiet. Adjustable tilt. 3 platform sizes. Stacking trays and multi-tiered versions save lab space.

NEXT>>>ADVANCE

BETTER WORKFLOW • ACCURATE RESULTS • HIGH VALUE

Next Advance, Inc. www.nextadvance.com

24 Prospect Avenue
Averill Park, NY 12018, USA
Telephone: 1.800.738.1681
Fax: 1.610.862.1718