

# THROW AWAY YOUR TIMERS!!!

## THE FREEDOM ROCKER™



*Model DX4 can process Westerns in up to 4 trays from blocking to final rinse.*

The **FREEDOM ROCKER™** automatically process blots and gels, including Westerns from blocker to final rinse. It is a rocker with a built-in pump and timer for transferring liquids to and from the trays. It frees you up from the tedious task of babysitting the blots or gels.

### AUTOMATED PROCESSING...

- > INCREASE PRODUCTIVITY
- > REDUCE DOWNTIME
- > SAVE MONEY

#### > INCREASE PRODUCTIVITY:

Make better use of your time. The **FREEDOM ROCKER™** automates the processing of blots and gels, freeing up time for more demanding activities. With accurate timing and pumping, results are more consistent.

#### > REDUCE DOWNTIME AND SHORTEN CYCLES:

THE **FREEDOM ROCKER™** allows you to process blots and gels anytime even during meetings and over weekends.

#### > SAVE MONEY: The **FREEDOM**

**ROCKER™** saves you money every time you use it, because you spend less money on paying someone to babysit a blot. Now they can do twice as much work for the same salary!

**RISK FREE: THE FREEDOM ROCKER™**  
comes with a 30 day money back guarantee

"I'm one of the first users of THE FREEDOM ROCKER™ at Amgen in Cambridge, MA. The machine gets LOTS of use from all of us here." — *Dr. Barbara Grubinska, Amgen I*

"It is perfect for silver staining of gels for mass spectrometry. It has made an UNBELIEVABLE impact on data generation for me in cancer protein screening." —*Dr. Steve Griffiths, Atlantic Cancer Research Institute*

"THE FREEDOM ROCKER™ is really convenient. I used it for washing gels while I worked on other things, and also as a regular rocker." —*Dr. Michele Lastro, The UAlbany Gen\*NY\*Sis Center for Excellence in Cancer*

## NEXT>>>ADVANCE

INNOVATIVE PRODUCTS • BETTER WORKFLOW  
ACCURATE RESULTS • [www.nextadvance.com](http://www.nextadvance.com)

24 Prospect Avenue  
Averill Park, NY 12018, USA  
Telephone: 800-738-1681  
Fax: 518-674-0189

# THROW AWAY YOUR TIMERS!!!

## AUTOMATICALLY PROCESS GELS/BLOTS.

The **FREEDOM ROCKER™ DX** automates the processing of gels and blots, using your same current protocols. Programming is extremely intuitive, with each program step the same as each protocol step. Program up to nine steps, where each step is an immersion in a reagent or a series of washes. For each step, simply set the volume and duration of the reagent or wash buffer, the number of dispenses, the rocking speed, and whether you want a clearing rinse. After the final wash, the buffer remains in the trays so that the gel or blot does not dry out

The **FREEDOM ROCKER™ DX** can be programmed for non-traditional protocols requiring initial rinses. It has the capacity to fill trays from 3 sources (one at a time to all trays), and to individually recover two reagents as well as aspirate into a waste container or sink.

Rocking speed up to 40 cycles/minute. Each programmable step is completely adjustable: time from 1min to 90hrs; volume from 5mL to 600mL; aspirate destination (to direct to waste bottle or reagent recovery bottle).

Reagent recovery gives you the opportunity to reuse precious antibodies, giving you the option for even more cost savings!

### Features of All FREEDOM ROCKERS™

- Automated delivery and aspiration of buffers/reagents.
- Reagent recovery for added cost savings.
- The FREEDOM ROCKER™ can call any lab bench home. Just plug into an electrical outlet.
- Quiet. No vacuum or air connections or vacuum pumps required.
- Drip proof quick-release connections.
- Patent pending technology.
- CE/UL certified 12V power supply (US or international).
- 2 year warranty.
- 30-day money back guarantee.

The **FREEDOM ROCKER DX4-MD** will deliver different reagents to each tray.



Model	Platform Size	Capacity	Special Features
<b>DX2</b>	10" x 12"	1 or 2 trays	
<b>DX2-MD</b>	10" x 12"	1 or 2 trays	multi-dispense -- different reagent for each tray
<b>DX4</b>	12" x 12"	4 trays; 1 or 2 w/ accessory kit	
<b>DX4-MD</b>	12" x 12"	4 trays; 1 or 2 w/ accessory kit	multi-dispense -- different reagent for each tray

**NEXT >>> ADVANCE**  
 INNOVATIVE PRODUCTS • BETTER WORKFLOW  
 ACCURATE RESULTS • [www.nextadvance.com](http://www.nextadvance.com)

24 Prospect Avenue  
 Averill Park, NY 12018, USA  
 Telephone: 800-738-1681  
 Fax: 518-674-0189