



Freedom Rocker Menu Flow Chart

Manual Run Settings Menu

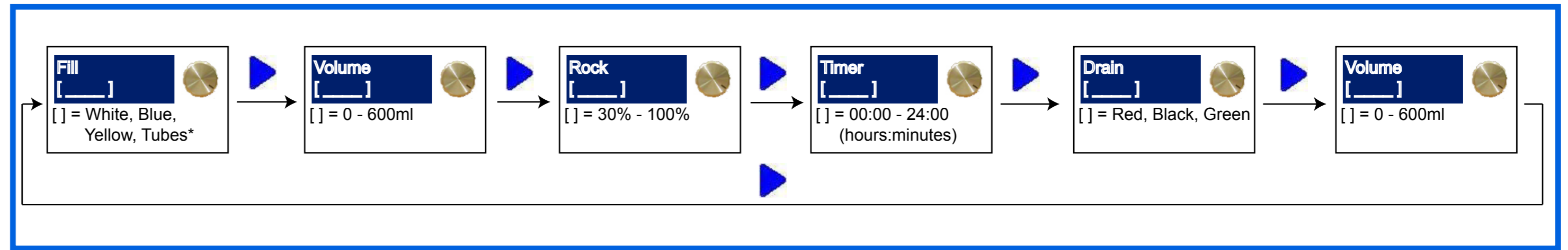


Automatic Run Settings Menu

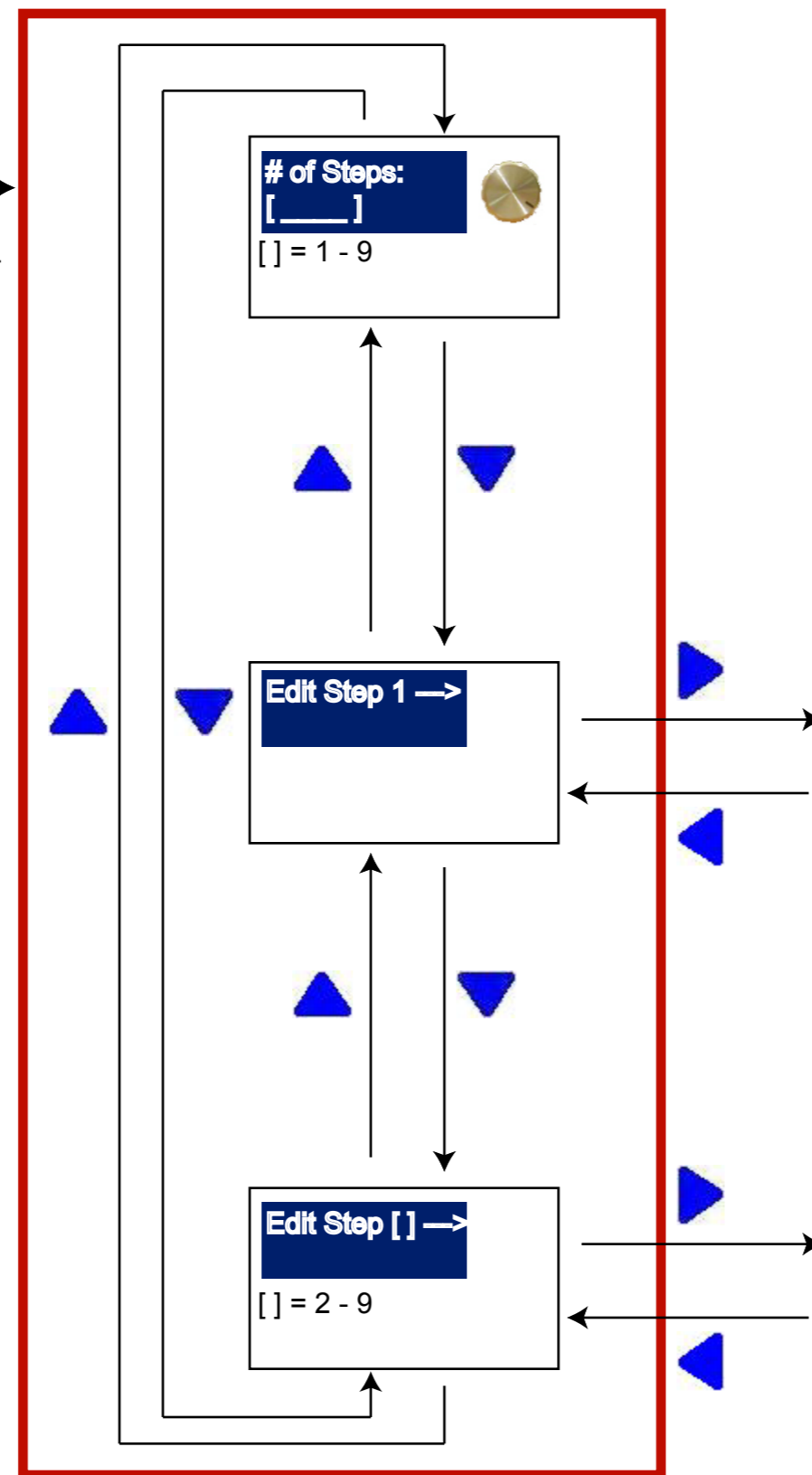

Notes

- The "Tubes" selection in the Fill menu only appears in multi-dispense models. This setting drains from the syringes on the dispense head.
- The two "volume" settings in the Manual Run Settings Menu both specify fill volume, and changing one will change both.
- At any point you may press "start" to begin the run. Pressing it again will pause the run. If you are in the Automatic Settings Menu or the Programming Menus, pressing start will run the automatic program. If you are in the Manual Run Settings Menu, pressing start will run the manual program.
- Pressing stop during the run will stop the run and return you to your previous position in the menu.
- Turning the knob during the run will change the rocking speed for that step

Program Step #1

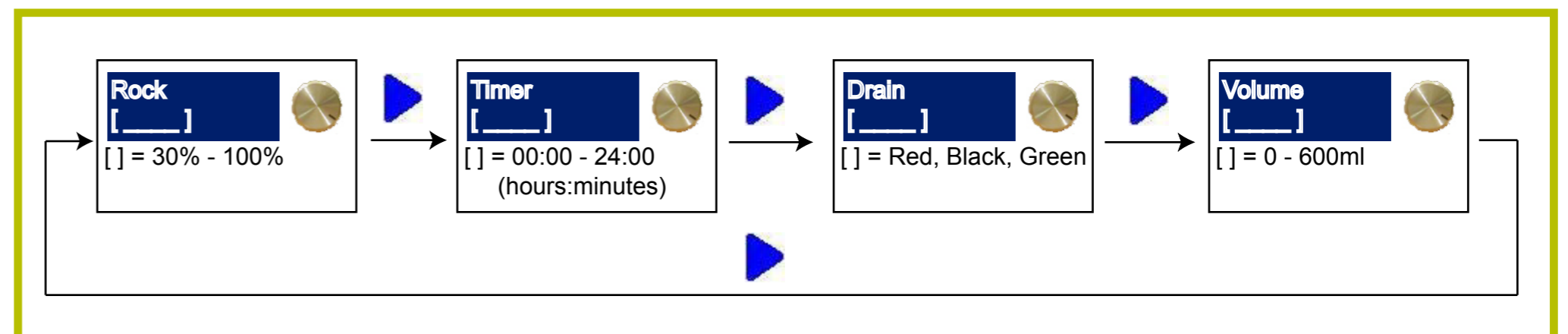

Program Step #2-9

