

# THROW AWAY YOUR TIMERS!!!

## THE FREEDOM ROCKER™



*Model DX4 can process Westerns in up to 4 trays from blocking to final rinse.*

The **FREEDOM ROCKER™** automatically processes Western, Northern, and Southern blots and gels from blocker to final rinse. It is a rocker with a built-in pump and timer for transferring liquids to and from the trays. It frees you up from the tedious task of babysitting the blots or gels.

- AUTOMATED PROCESSING...**
- > INCREASE PRODUCTIVITY
  - > REDUCE DOWNTIME
  - > SAVE MONEY

**> INCREASE PRODUCTIVITY:**

Make better use of your time. The **FREEDOM ROCKER™** automates the processing of blots, gels and slides, freeing up time for more demanding activities. With accurate timing and pumping, results are more consistent.

**> REDUCE DOWNTIME AND SHORTEN CYCLES:**

The **FREEDOM ROCKER™** allows you to process blots, gels and slides anytime - even during meetings and over weekends.

**> SAVE MONEY:** The **FREEDOM**

**ROCKER™** saves you money every time you use it, because no scientist is required to babysit the blot/gel. Now they can do twice as much work for the same salary!

***RISK FREE: THE FREEDOM ROCKER™ comes with a 30 day money back guarantee***

"I'm one of the first users of THE FREEDOM ROCKER™ at Amgen in Cambridge MA. The machine gets LOTS of use from all of us here." – *Dr. Barbara Grubinska, Amgen*

"It is perfect for silver staining of gels for mass spectrometry. It has made an UNBELIEVABLE impact on data generation for me in cancer protein screening." – *Dr. Steve Griffiths, Atlantic Cancer Research Institute*

"THE FREEDOM ROCKER™ is really convenient. I used it for washing gels while I worked on other things, and also as a regular rocker." – *Dr. Michele Lastro, The UAlbany Gen\*NY\*Sis Center for Excellence in Cancer Genomics*

**NEXT >>> ADVANCE**

INNOVATIVE PRODUCTS • BETTER WORKFLOW  
ACCURATE RESULTS • [www.nextadvance.com](http://www.nextadvance.com)

24 Prospect Avenue  
Averill Park, NY 12018, USA  
Telephone: 800-738-1681  
Fax: 518-674-0189  
[info@nextadvance.com](mailto:info@nextadvance.com)

# THROW AWAY YOUR TIMERS!!!

## AUTOMATICALLY PROCESS GELS/BLOTS/SLIDES

Are you chained to your timer? Throw it out and use the **FREEDOM ROCKER™** to automate your work. Simply place your blots, gels or slides in trays, set the program, connect the bottles/tubes of reagents and walk away. The microcontroller ensures precise and consistent results. Multi-dispense models can even dispense different reagents to each tray.

The **FREEDOM ROCKER™** can run in manual mode to perform a dispense, timed rocking, or aspiration step, or in automatic mode where it can perform up to 9 protocol steps. For each step, simply set the volume of the reagent or wash buffer, the duration, the number of dispenses, the rocking speed, and whether you want a clearing rinse. After the final wash, the buffer remains in the trays so that the gel or blot does not dry out. For each step, select among 3 buffers/reagents and aspirate to waste container or either of 2 recovery containers.

Rocking speeds up to 40 cycles/minute. Each programmable step is completely adjustable: time from 1 min to 90hrs; volume from 5mL to 600mL; aspirate destination direct to waste bottle or reagent recovery bottle.

Reagent recovery gives you the opportunity to reuse precious antibodies for even more cost savings!

### Features of All **FREEDOM ROCKERS™**

- Automated delivery and aspiration of buffers/reagents.
- Reagent recovery for added cost savings.
- The **FREEDOM ROCKER™** can call any lab bench home. Just plug into an electrical outlet.
- Quiet- No air connections or vacuum pumps required.
- Drip proof quick-release connections.
- Patent pending technology.
- CE/UL certified 12V power supply (US or international).
- 2 year warranty.
- 30-day money back guarantee.

The **FREEDOM ROCKER™ DX4-MD** will deliver different reagents to each tray.



Model	Platform Size	Capacity	Special Features
<b>DX2</b>	10" x 12"	1 or 2 trays	
<b>DX2-MD</b>	10" x 12"	1 or 2 trays	multi-dispense -- different reagent for each tray
<b>DX4</b>	12" x 12"	4 trays; 1 or 2 w/ accessory kit	
<b>DX4-MD</b>	12" x 12"	4 trays; 1 or 2 w/ accessory kit	multi-dispense -- different reagent for each tray

## NEXT >>> ADVANCE

INNOVATIVE PRODUCTS • BETTER WORKFLOW  
ACCURATE RESULTS • [www.nextadvance.com](http://www.nextadvance.com)

24 Prospect Avenue  
Averill Park, NY 12018, USA  
Telephone: 800-738-1681  
Fax: 518-674-0189  
[info@nextadvance.com](mailto:info@nextadvance.com)