

Operator's Manual




Next Advance Rockers

Congratulations!

Congratulations on your purchase of a Rocker from Next Advance. Please read this operator's manual which explains proper operation of the instrument. This manual can be viewed on the web at https://www.nextadvance.com/pdf/Rocker_Manual.pdf.


You can view our other products at www.nextadvance.com.

PARTS OF THE ROCKER

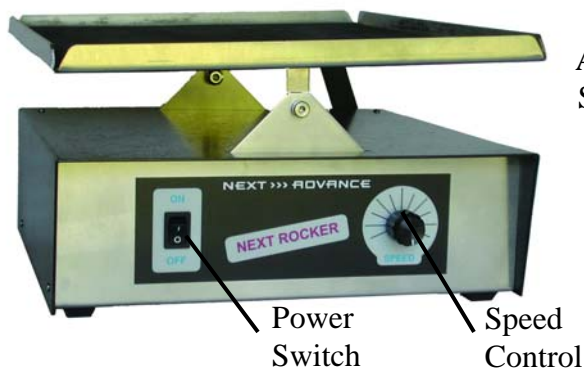

Front of Rocker

Tilt
Adjust
Screw

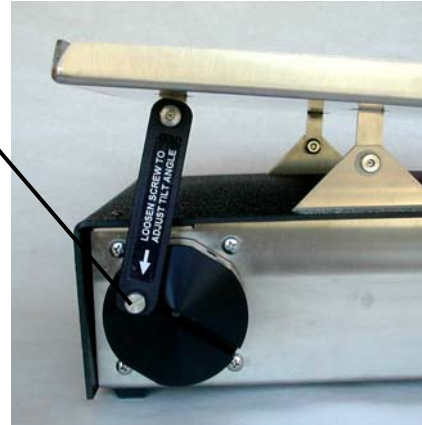

Rear of Rocker

SETUP

Place the Rocker on a level lab bench. Plug the power supply into the back of the Rocker and then plug it into a wall outlet. It is now set up.

OPERATION

Turn on the Power Switch and turn the Speed Control knob to adjust the rocking rate.

To adjust the tilt angle, use a flat blade screwdriver to loosen the labeled Tilt Adjust Screw on the rear of the instrument, slide the linkage to the desired position, then tighten the screw.

TROUBLESHOOTING

In addition to the tips given below, a thorough list of troubleshooting tips is at <https://www.nextadvance.com/content/infinity-rocker-next-rocker-troubleshooting-116.cfm>.

If nothing happens, the power supply plug might not be in a live wall outlet or the power supply connector plug might not be fully in the jack on the back of the Rocker.

If the unit stops working, turn the system off for 15 minutes to allow the electronics to reset. If the Rocker does not turn on after this period, contact customer service.

SUPPORT

A thorough list of FAQs, is at <https://www.nextadvance.com/product/rocker-laboratory-platform-shaker-214.cfm/#tab3/top>.

If you cannot find an answer there, please contact customer service by email at support@nextadvance.com or by telephone.

WARRANTY

Next Advance warrants its Products against defects in materials and workmanship for time periods which vary according to the Product. Within these time periods, Next Advance will replace or repair, without charge to the original purchaser, any part which is defective.

Next Rocker	One year
Infinity Rocker	Four years

The warranty is void if the Product is defective due to product accident, product modification, exposure to radiation other than for sterilization, connection to an improper electrical supply, lack of proper maintenance, contamination, improper installation or misuse. The warranty shall also not apply to defects arising from fire, flood, lightning or other conditions unrelated to correct operation of the Product.

OPERATOR'S MANUAL FOR ROCKERS

Next Advance's liability is limited, at the company's election, to (1) refund of the original purchaser's purchase price for the Product (2) repair of the Product, or (3) replacement of the Product or defective parts. Evidence of purchase by the original purchaser is required. Next Advance may also request documentation of proper maintenance, if applicable.

Next Advance makes no other warranty, express or implied, with respect to its Products. NEXT ADVANCE MAKES NO WARRANTY RESPECTING THE MERCHANTABILITY OF THE PRODUCTS OR THEIR SUITABILITY OR FITNESS FOR ANY PARTICULAR PURPOSE OR USE. Next Advance shall not be liable for, indirect, special, incidental or consequential damages of any nature. Any recovery for any claim shall be limited to the original purchase price for the product.

Operator's Responsibility

Provide proof of purchase and provide normal care and maintenance.